

# Spin-X® UF Concentrators FAQs

CORNING



1. Can I re-use a Spin-X UF concentrator?
2. Can I use the Spin-X UF devices with a swinging-bucket rotor?
3. Can I use the molecular weight cut-off (MWCO) of Spin-X UF devices as a guide for concentrating DNA?
4. Can the Spin-X UF concentrators be spun to dryness?
5. Do I need to pre-rinse the Spin-X UF concentrators before I add my sample?
6. How can I increase my recovery with the Spin-X UF concentrators?
7. How can I tell what the MWCO of my Spin-X UF concentrator is?
8. How long does it take to concentrate a sample using Spin-X UF concentrators?
9. How much of the protein in my sample can I expect to recover with the Spin-X UF concentrators?
10. Is it possible to concentrate and desalt my protein sample using the Spin-X UF concentrators?
11. What rotor sizes are compatible with the Spin-X UF concentrators?
12. What is the housing material of the Spin-X UF concentrators?
13. Do you have any recommendations in which direction the Spin-X UF device should be placed in the centrifuge rotor?
14. What membranes are available in the Spin-X UF concentrators?
15. What pipet or pipet tips do you recommend for recovering the concentrate from the Spin-X UF concentrator?
16. Can Spin-X UF concentrators be sterilized with ethanol?
17. Can I autoclave the Spin-X UF concentrators?
18. What is the purpose of the small hole centered in the UF 6 and 20 caps?
19. What is the maximum sample volume for the different Spin-X UF concentrators?
20. What is the maximum speed to spin the different Spin-X UF concentrators?
21. What is the membrane surface area in the Spin-X UF concentrators?
22. Can Spin-X UF concentrators be used for fractionation?

## Spin-X® UF Concentrators Frequently Asked Questions

### 1. Can I re-use a Spin-X UF concentrator?

No, the device is developed for one time use only.

### 2. Can I use the Spin-X UF devices with a swinging-bucket rotor?

All Spin-X UF (except Spin-X UF 500) can be used in swinging-bucket rotors. The vertical membrane gives fast flow rates and the sample can not be centrifuged to dryness, due to the integrated dead stop volume.

### 3. Can I use the molecular weight cut-off (MWCO) of Spin-X UF devices as a guide for concentrating DNA?

No, concentrating DNA, which has a linear structure, requires that Spin-X UF membranes be characterized in terms of a base pair cut-off as indicated by the following table.

MWCO	Ds DNA base pairs
5,000	15
10,000	30
30,000	70
50,000	150
100,000	200

### 4. Can the Spin-X UF concentrators be spun to dryness?

No, because all Spin-X UF concentrators have an integrated dead stop volume and there is no risk to spin the sample to dryness.

### 5. Do I need to pre-rinse the Spin-X UF concentrators before I add my sample?

Membranes in the Spin-X UF concentrators contain trace amounts of glycerine and sodium azide. These compounds can be removed by rinsing the device with buffer solution or deionised water. Decant filtrate and concentrate before processing sample solution. If a device is to be prepared for later use, make sure the membrane does not dry out by storing the device at 4°C with deionized water or buffer.

### 6. How can I increase my recovery with the Spin-X UF concentrators?

If you have a low concentration protein sample, we recommend passivating (blocking) the membrane surface. The intention of the passivation procedure is to improve recovery of protein samples by pretreating the device. For this purpose a range of solutions are suggested in the the next table. For more detailed information, refer to the application note.

Type	Solution
Powdered milk	1% in ultra pure water
BSA	1% in PBS
Tween 20	5% in ultra pure water
SDS	5% in ultra pure water
Triton-X 100	5% in ultra pure water
PEG 3000	5% in ultra pure water

### 7. How can I tell what the MWCO of my Spin-X UF concentrator is?

This information is printed on the device.

### 8. How long does it take to concentrate a sample using Spin-X UF concentrators?

It depends on the starting conditions, but for example to concentrate 1mg BSA 30 x at 20°C in a swing-bucket rotor with a 30,000 MWCO device:

Device	Minutes
Spin-X UF 500	5
Spin-X UF 6	12
Spin-X UF 20	22

### 9. How much of the protein in my sample can I expect to recover with the Spin-X UF concentrators?

If you choose the correct MWCO, >90% recovery.

### 10. Is it possible to concentrate and desalt my protein sample using the Spin-X UF concentrators?

Yes, it is possible. First you concentrate your protein; then you refill the device with the buffer containing a lower salt concentration and spin again. After this step, you have desalted and concentrated your protein. For more information please refer to our desalting application note.

**11. What rotor sizes are compatible with the Spin-X UF® concentrators?**

- Spin-X UF 500 can be processed in all fixed angle rotors accepting 2.2 mL tubes.
- Spin-X UF 6 can be processed in all rotors accepting 15 mL conical bottom tubes.
- Spin-X UF 20 can be processed in all rotors accepting 50 mL conical bottom tubes.

**12. What is the housing material of the Spin-X UF concentrators?**

The housing material is polycarbonate.

**13. Do you have any recommendations in which direction the Spin-X UF device should be placed in the centrifuge rotor?**

- If you use a fixed angle rotor, the Spin-X UF concentrator must be placed so that the graduation faces outwards/upwards.
- If you use a swing bucket rotor, you do not need to place the Spin-X UF concentrator into the rotor in any specific way.

**14. What membranes are available in the Spin-X UF concentrators?**

All Spin-X UF concentrators have a PES (Polyethersulfone) membrane.

**15. What pipet or pipet tips to you recommend for recovering the concentrate from the Spin-X UF concentrator?**

We recommend an adjustable pipettor with a yellow tip or gel loader tips for optimal recovery.

**16. Can Spin-X UF concentrators be sterilized with ethanol?**

Yes, use a 70% ethanol solution or sterilising gas mixture.

**17. Can I autoclave the Spin-X UF concentrators?**

The Spin-X UF concentrators cannot be autoclaved.

**18. What is the purpose of the small hole centered in the UF 6 and 20 cap?**

The small hole in the UF 6 and 20 caps serves to equalize pressure differentials during centrifugation.

**19. What is the maximum sample volume for the different Spin-X UF concentrators?**

Device	Max. volume
Spin-X UF 500	500 µL
Spin-X UF 6	6 mL
Spin-X UF 20	20 mL

**20. What is the maximum speed to spin the different Spin-X UF concentrators?**

Concentrator	Maximum Recommended Centrifugal Force		
	Spin-X UF 500	Spin-X UF 6	Spin-X UF 20
<i>Maximum Spin Force – Swing Bucket</i>			
5,000 to 50,000 MWCO PES	Do not use	4,000 x g	5,000 x g
>100,000 MWCO PES	Do not use	4,000 x g	3,000 x g
<i>Maximum Spin Force – Fixed Angle</i>			
5,000 to 50,000 MWCO PES	15,000 x g	10,000 x g	8,000 x g
>100,000 MWCO PES	15,000 x g	6,000 x g	6,000 x g

**21. What is the membrane surface area in the Spin-X UF concentrators?**

Device	Active membrane surface area
Spin-X UF 500	0.5 cm <sup>2</sup>
Spin-X UF 6	2.5 cm <sup>2</sup>
Spin-X UF 20	6.0 cm <sup>2</sup>

**22. Can Spin-X UF concentrators be used for fractionation?**

Ultrafiltration cannot be used for fractionation. The ultrafiltration membranes do not allow separation of proteins from a protein mixture, as proteins smaller than the cut-off will not necessarily pass the membrane completely.

## Corning® Spin-X® UF Concentrator Ordering Information

Cat. No.	Description	Capacity	Membrane	Pack Size
431477	Spin-X UF 500	500 µL	5,000 MWCO	25
431478	Spin-X UF 500	500 µL	10,000 MWCO	25
431479	Spin-X UF 500	500 µL	30,000 MWCO	25
431480	Spin-X UF 500	500 µL	50,000 MWCO	25
431481	Spin-X UF 500	500 µL	100,000 MWCO	25
431482	Spin-X UF 6	6 mL	5,000 MWCO	25
431483	Spin-X UF 6	6 mL	10,000 MWCO	25
431484	Spin-X UF 6	6 mL	30,000 MWCO	25
431485	Spin-X UF 6	6 mL	50,000 MWCO	25
431486	Spin-X UF 6	6 mL	100,000 MWCO	25
431487	Spin-X UF 20	20 mL	5,000 MWCO	12
431488	Spin-X UF 20	20 mL	10,000 MWCO	12
431489	Spin-X UF 20	20 mL	30,000 MWCO	12
431490	Spin-X UF 20	20 mL	50,000 MWCO	12
431491	Spin-X UF 20	20 mL	100,000 MWCO	12

For additional product or technical information, please visit [www.corning.com/lifesciences](http://www.corning.com/lifesciences) or call 1.800.492.1110. Outside the United States, please call +1.978.442.2200.

# CORNING

### Corning Incorporated Life Sciences

Tower 2, 4th Floor  
900 Chelmsford St.  
Lowell, MA 01851  
t 800.492.1110  
t 978.442.2200  
f 978.442.2476

[www.corning.com/lifesciences](http://www.corning.com/lifesciences)

### Worldwide Support Offices

#### ASIA / PACIFIC

**Australia/New Zealand**  
t 65 6733-6511  
f 65 6735-2913

**China**  
t 86-21-5467-4666  
f 86-21-5407-5899

**India**  
t 91-124-235 7850  
f 91-124-401 0207

#### Japan

t 81 (0) 3-3586 1996/1997  
f 81 (0) 3-3586 1291/1292

#### Korea

t 82 2-796-9500  
f 82 2-796-9300

#### Singapore

t 65 6733-6511  
f 65 6861-2913

#### Taiwan

t 886 2-2716-0338  
f 886 2-2716-0339

#### EUROPE

##### France

t 0800 916 882  
f 0800 918 636

##### Germany

t 0800 101 1153  
f 0800 101 2427

##### The Netherlands

t 31 20 655 79 28  
f 31 20 659 76 73

##### United Kingdom

t 0800 376 8660  
f 0800 279 1117

### All Other European Countries

t 31 (0) 20 659 60 51  
f 31 (0) 20 659 76 73

#### LATIN AMERICA

##### Brasil

t (55-11) 3089-7419  
f (55-11) 3167-0700

##### Mexico

t (52-81) 8158-8400  
f (52-81) 8313-8589